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Cat: G1355

Size: 3×50mL

Storage: RT, avoid light, valid for 1 year.

Kit Components

Reagent	3×50mL	Storage
Reagent A: Mallory Stain A	50mL	RT , avoid light
Reagent B: Mallory Stain B	50mL	RT , avoid light
Reagent C: Mallory Stain C	50mL	RT , avoid light

Introduction

Trichrome Stain usually refers to the staining of nucleus and the selective stain of collagen and muscle fibers. The dyeing principle of this method is related to the size of anionic dye molecules and the permeability of the tissues: the size of the molecules is reflected by the molecular weight, and the small molecular weight can easily penetrate the compact and low permeability tissues, while the large molecular weight can only enter the loose and high permeability tissues.

Mallory Trichrome Stain Kit can dye collagen fiber blue, nucleus and muscle fiber red, and red blood cell orange.

Protocol(for reference only)

1. Dewax paraffin sections and rehydrate in graded of alcohol.
2. Stain with Mallory Stain A for 1-3mins.
3. Stain with Mallory Stain B for 2mins.
4. Stain with Mallory Stain C for 15mins.
5. Differentiate by 95% alcohol for 10-15s (control under microscope).
6. Conventional dehydration, transparentize by xylene.
7. Seal with resinene.

Result

Nucleus	Red
Red blood cell	Orange
Muscle fiber	Red
Collagen fiber	Blue

Note

1. In the dyeing process, go directly to the next step after 2/3/4 steps without washing.
2. For your safety and health, please wear experimental clothes and disposable gloves.