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规格: 100mL/500mL

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Cat: G1450

Size: 100mL/500mL

Storage: 2-8°C, avoid light, valid for 6 months.

Introduction

Calcium is abundant in human body. As a scaffold to support human body, bone plays an important role in secretion, transportation, muscle contraction, nerve conduction and so on. Alizarin Red S is an anthraquinone derivative, which is sodium alizarin sulfonate. It can chelate with calcium carbonate or calcium phosphate to form orange red complex. Generally speaking, the effect of these dyes is better than that of slight staining when they are used to recognize large amount of calcium. But Alizarin Red S can get more reliable results for a small amount of sediment. It is often combined with fast green or Mayer hematoxylin staining solution to form orange red precipitate, which is suitable for dyeing a small amount of calcium salt tissue.

Protocol(for reference only)

For paraffin section

1. Fix the tissue in 10% neutral formalin or ethanol, then dehydrate and embed.
2. Place the section into 95% ethanol.
3. Place the section vertically and air dry thoroughly.
4. Add the section into a vat containing Alizarin Red S Solution, 0.2%, pH8.3 and stain for 5-10min.
5. Quickly rinse in distilled water.
6. (optional)Re-dyeing with counterstain solution and wash with distilled water for three times.
7. Conventionally dehydrate, transparent and seal with resinene.

For cultured cell

1. Remove the medium from the plate and wash twice with PBS.
2. Fix in 10% neutral formalin or 4% paraformaldehyde for 10-15min.
3. Discard the fixative and wash with distilled water for 3 times.
4. After the water is completely absorbed, add Alizarin Red S Solution, 0.2% , pH8.3 slowly and dye for 20-30min.
5. Discard the dye and wash 3-5 times with distilled water.
6. Add a proper amount of distilled water into each hole to prevent drying in the hole. Observe and photograph under the microscope.

Result

Calcium Deposits	Orange Red
------------------	------------

Note

1. The dyeing time of Alizarin Red S Solution should be determined according to the content of calcium salt. It should be observed under the microscope. When the calcium salt is deep orange red, take out and wash.
2. After dyeing with Alizarin Red S Solution, the calcium deposits are birefringent.
3. When re-dyeing with fast green, the background is green. When re-dyeing with Mayer's hematoxylin staining solution, the nucleus is blue.
4. This method is particularly useful in identifying and detecting small amounts of calcium, such as abnormal calcium in the kidney (excessive calcium in the urine).
5. For your safety and health, please wear experimental clothes and disposable gloves.